










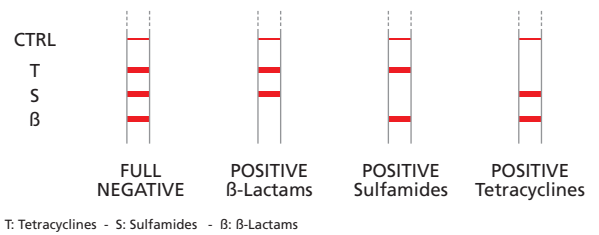
TRISENSOR is a Rapid Test for simultaneous detection of β -lactams, Tetracyclines and Sulfonamides in milk in 6 minutes.

Benefits


- 
 → **MULTIPLE**
 Detection of β -lactams, Tetracyclines and Sulfonamides.
- 
 → **RAPID**
 Results in 6 minutes.
- 
 → **SENSITIVE**
 EU MRLs.
- 
 → **RELIABLE AND ROBUST**
- 
 → **USER-FRIENDLY**
 Easily performed on site or in lab.
- 
 → **COST-EFFECTIVE**
- 
 → **ECOLOGICAL PACKAGING**

Results

Visual interpretation of the result by comparing the intensity of the "TEST" lines with the one of the "CTRL" line.



How to use

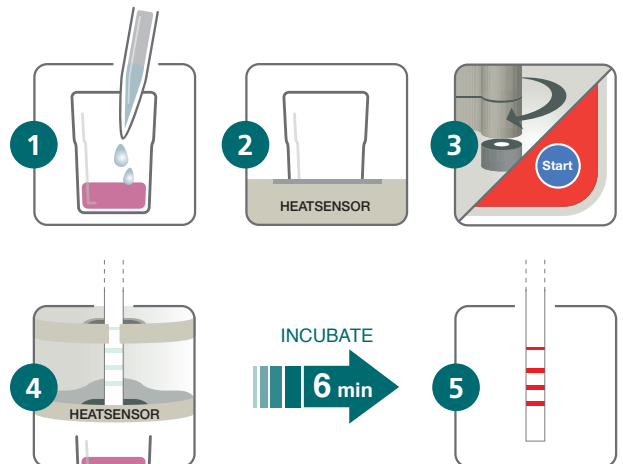


→ 3 CLASSES AT A TIME

→ β -LACTAMS + TETRACYCLINES + SULFONAMIDES

Sensitivity

β -LACTAMS	LOD (PPB)	SULFONAMIDES	LOD (PPB)	TETRACYCLINES	LOD (PPB)
PENICILLINS		Sulfadiazine	5-6	Tetracycline	80-100
Penicillin-G	2-3	Sulfapyridine	0,5-1	Oxytetracycline	50-60
Ampicillin	3-5	Sulfathiazole	4-8	Chlortetracycline	50-55
Amoxicillin	3-5	Sulfamethoxazole	150-300	Doxycycline	20-30
Oxacillin	12-18	Sulfamethazine	1-1,5		
Cloxacillin	6-8	Sulfamethoxypyridazine	1-3		
Dicloxacillin	6-8	Sulfadimethoxine	10-15		
Nafcillin	30-40	Sulfacetamide	300-600		
CEFALOSPORINS		Sulfamerazine	1,5-2,5		
Ceftiofur	10-15	Sulfamonomethoxine	7-14		
Cefquinome	30-35	Sulfaquinoxaline	15-30		
Cefazolin	18-22	Sulfachloropyridazine	5-10		
Cefapirine	6-8	Sulfaguanidine	15-25		
Cefacetrile	30-40	Sulfamethizole	75-100		
Cefoperazone	3-4	Sulfasalazine	250-350		
Cefalexin	1000-1200				
Cefalonium	3-5				

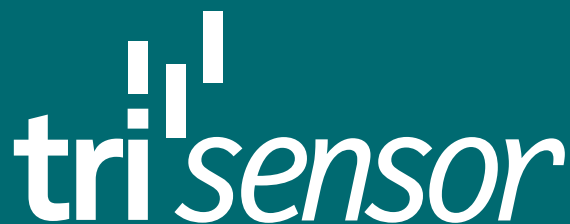

Products References

KIT035 → TRISENSOR MK BTS MRL - 96 pc

APP032 → HEATSENSOR DUO

APP039 → READSENSOR ESE FULL CASE

unisensor 

Unisensor s.a.

ZI du Dossay
Rue du Dossay, 3
B-4020 Wandre (Liège)
BELGIUM

Phone +32 4 252 66 02
Fax +32 4 252 90 55
info@unisensor.be

www.unisensor.be

Reseller

NOACK
GROUP OF COMPANIES

Sales & Service: www.noackgroup.com